



EI COULBOURN INSTRUMENTS

CLEANING ACT-55X SERIES CAGES

The cages used for the ACT-551, ACT-551MS, ACT-552, and ACT-552MS can be autoclaved. Allentown Caging recommends using chlorine dioxide however the cages must be rinsed with clean water. If autoclaved with chlorine dioxide still on the cage, residue may stick to the cage and turn white. If the cages do not need to be sterilized, the cages can be washed with a regular detergent. Be sure the switch assembly is removed from the cage before washing (these just pull out and can be pushed back in). There is a lock washer used on the cage assembly that has a temperature limit of 250 degrees F so if autoclaved, the temperature must not exceed 250 degrees F.